

# Quick Installation Guide for Pall Allegro™ 3D Biocontainers – 100, 200, 500, 1000 and 1500 L



1

■ If trolley or casters are being used, ensure brakes are on



2

■ If using levelling feet, adjust levels while empty



3

■ Open the doors of the tote




4

■ Open package. Confirm Allegro system is not damaged and locating plate is in place on the bottom port




5

■ Confirm tubing on inlet and outlet ports is securely connected

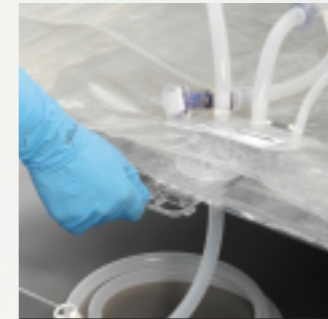

6

■ Slide the Allegro biocontainer into the tote and align the locating plate in the cutaway in the bottom of the tote

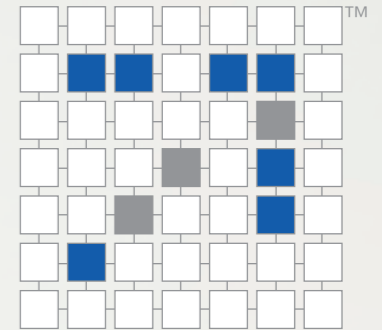


**AllegroSystems**  
The Single-Use Solution



7

■ Ensure side panels of the Allegro biocontainer are facing upwards, with tubing and components connected to the top ports placed in the center of the folded biocontainers

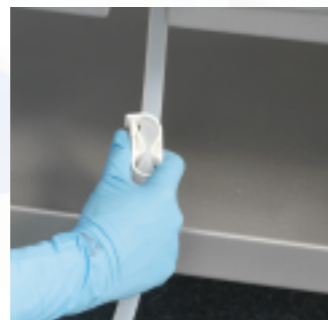

8

■ Check all outlet lines are closed  
■ Open filling lines



9

■ Close and lock the door



10

■ Start filling



11

■ Once filling is complete, close the filling line  
■ For dispensing, open the outlet line. Fluid can be dispensed by gravity or with the use of a pump



12

■ Once Allegro biocontainer is empty, discard the Allegro system in accordance with instructions at your facility



■ Stack 200 L totes (maximum 3) and 500, 1000 and 1500 L totes (maximum 2)



■ **DO NOT** lift totes manually



■ **DO NOT** stack the totes on a trolley or casters



■ **DO NOT** forcibly unfold Allegro biocontainer prior to use



■ **DO NOT** open the door while Allegro biocontainer is filling or full



■ **DO NOT** overfill Allegro biocontainers over their maximum volume specification