


Quick Installation Guide

Kleenpak™ Sterile Disconnectors

Component Parts

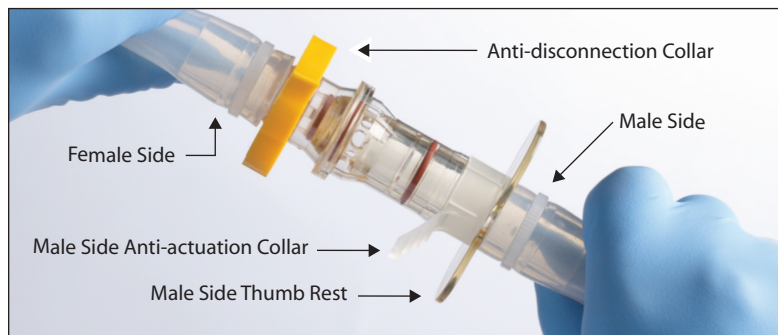

Four Simple Steps to Operate



Step 1: Drain as much fluid as possible from the disconnector, clamp off tubing on both sides of the device and pull off (remove) male side anti-actuation collar.



Step 2: Push male section fully into female section until audible 'click' and male side cannot be pushed any further.

⚠ See Warning



Step 3: Pull off (remove) anti-disconnection collar towards tubing.

Additional Information:

For more detailed instructions on how to operate the Kleenpak Sterile Disconnector, please refer to Pall publication USD2727, which is included in each box of disconnectors and is available from Pall on request.

⚠ Warning: Activation must take place when the connector is either free of liquid or the female side tubing is open and unrestricted.



Step 4: Twist, then separate the two sections.



Optional Step:
Fit female end cover to disconnected female section.

Corporate Headquarters

Port Washington, NY, USA
+1 800 717 7255 toll free (USA)
+1 516 484 5400 phone
biopharm@pall.com e-mail

European Headquarters

Fribourg, Switzerland
+41 (0)26 350 53 00 phone
LifeSciences.EU@pall.com e-mail

Asia-Pacific Headquarters

Singapore
+65 6389 6500 phone
sgcustomerservice@pall.com e-mail

Visit us on the Web at www.pall.com/allegro

E-mail us at allegro@pall.com

International Offices

Pall Corporation has offices and plants throughout the world in locations such as: Argentina, Australia, Austria, Belgium, Brazil, Canada, China, France, Germany, India, Indonesia, Ireland, Italy, Japan, Korea, Malaysia, Mexico, the Netherlands, New Zealand, Norway, Poland, Puerto Rico, Russia, Singapore, South Africa, Spain, Sweden, Switzerland, Taiwan, Thailand, the United Kingdom, the United States, and Venezuela. Distributors in all major industrial areas of the world. To locate the Pall office or distributor nearest you, visit www.pall.com/contact.

The information provided in this literature was reviewed for accuracy at the time of publication. Product data may be subject to change without notice. For current information consult your local Pall distributor or contact Pall directly.