Pall® Kleenpak™ Water Filter

ENGLISH

KLEENPAK™ WATER FILTER

Re-order No. KAQ14F1S, KAQ14F1R

DESCRIPTION
The Kleenpak Water Filter is supplied sterilized and integrity tested. The sterilizing gase "Supor" membrane within the filter cartridge is composed of a 0.2 µm polypropylene membrane for the removal of bacteria, protozoa, fungi and particles* from water used for hand hygiene and particles* present in a food outlet, as well as for the washing of hands in the saas des salles blanches des industries pharmaceutiques ainsi que pour les applications similaires. Il est destiné à un usage unique.

STERILE
Sterilized by gamma irradiation.

INSTALLATION

(Read and follow carefully)

Recommendations for filter installation, positioning and use must be followed to minimize inadvertent microbial contamination downstream of the filter membrane.

1. Filter Attachment to Water Outlet Point

Ensure O-ring in position to groove before attaching filter to water outlet point. The filter should be fitted using a quick connect adaptor, which allows rapid attachment and removal of the filter to taps or to the end of a shower hose**.

The filter may be used on a water outlet point to provide water for applications described above.

The filter should be positioned to avoid or minimize back splash to the outlet when in use.

The Kleenpak Water Filter should be handled carefully during installation and touching the filter outlet during use should be avoided.

The water should be positioned directly into the sink drainage point.

The filter outlet should not be immersed in collected water.

The Kleenpak Water Filter contains an effective bacteriostatic compound to minimize the risk of inadvertent retrograde contamination of the filtrant.

Depend on local environmental and hygiene factors.

The external joints of the filter and connectors may be cleaned using sanitizing agents containing iso-propyl alcohol, quaternary ammonium compounds, benzyl ammonium chlorides, iodides and non-ionic surfactants. Periodic wiping or spraying of these sanitizing agents onto the filter housing or housing outlet does not affect filter integrity. At no point should the filter be disconnected during this procedure.

Always routinely water system disinfection the filter can be left in place to enable the outlet to be flushed but a new filter must be installed immediately after flushing and before the outlet is brought back into service.

Reduction of water flow may occur as a result of heavy contamination of the water with particles, native micro-organisms. This may reduce the filter's lifetime. If this occurs, please inform your local Pall representative or distributor, who may advise on suitable pre-filtration.

PRECAUTIONS

The filter is designed for use at a maximum upstream operating pressure of 5 bar (approx. 75 psi), at a maximum influent water temperature of 60 °C. The filter membrane can withstand a maximum continuous operating pressure of 10 bar (150 psi) for a period of 30 minutes over the life of the filter. Whilst in use, the filter should not be used to supply water to the downstream side and/or entrapment into the filter which may cause air-locking and significant reduce water flow. If leakage is observed connection to the water outlet, the filter should be removed and replaced with a new filter. The cause of the leakage should be investigated and corrective action taken. Ensure that fitting of the Kleenpak Water Filter is compliant with local water regulations. Avoid mechanical impact as this may damage the filter and/or connector. Used filters should be disposed of according to local procedures.

** See product literature for performance summary.

** For information on suitable adaptors contact your local Pall representative.

PALL, Pall, Kleenpak and Supor are trade marks of Pall Corporation. Filtration, Separation, Solution is a service mark of Pall Corporation.

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PALL Pall, Kleenaak ja Supor on firma Pall Corporation kaubamärgid.
Filtration, Separation, Solution - a Pall Corporation business.