



## Protocol for Genomic DNA Purification from Fresh Cultured CHO (Chinese Hamster Ovary) Cells Using Pall Nucleic Acid Binding Nanosep® Centrifugal Device

### 1. Consumables and Reagents

**Table 1**

Consumables for gDNA Purification (nuclease-free consumables are recommended)

Supplier	Product Description	Part Number
Pall Laboratory	Nucleic Acid Binding (NAB) Nanosep Centrifugal Device	ODNABC33, ODNABC34
VWR	Ethanol (not denatured)	71001-866
VWR	Spectrophotometer Cuvettes ~100 µL (260/280 nm)	47743-840
VWR	Tubes 15 mL (RNase-DNase free)	89401-574
VWR	Tubes 50 mL (RNase-DNase free)	89401-572
VWR	Microcentrifuge Tubes 1.5 mL (RNase-DNase free)	76005-210

**Table 2**

Reagents for gDNA Purification (nuclease-free reagents are recommended)

Supplier	Product Description	Part Number
Qiagen	Buffer AL (264 mL)	19075
Qiagen	Buffer ATL (200 mL)	19076
Qiagen	Buffer AW1 (242 mL)	19081
Qiagen	Buffer AW2 (324 mL)	19072
Qiagen	Proteinase K	19131
Qiagen	RNase A (100 mg/mL)	19101
Qiagen	Nuclease-free water	129115
VWR	Tris Buffer pH 7.0 (1 M)	89500-584
VWR	PBS Tablet	97062-732

### 2. Instruments

- Microcentrifuge
- Spectrophotometer
- Vortex
- Water bath or heating block

### 3. Important Points Before Starting

- Buffer AW1 and Buffer AW2 are supplied as concentrates. Before using for the first time, add the appropriate amount of ethanol (96-100 %) as indicated on the bottle to obtain a working solution. Before each use, mix Buffer AW1 by inverting several times.
- All centrifugation steps are performed at room temperature at 10,000 – 14,000 x g.
- Use only the collection tubes provided in the NAB Nanosep device box. There is enough to complete this process in full.
- All buffers should be allowed to equilibrate to room temperature before use.
- Briefly centrifugate the tubes after vortexing to remove drops from inside the lid.
- Increase volumes of lysis buffers and reagents proportionally for samples >200 µL.
- Ensure that any precipitate formed during any of the below steps is not lost.
- Centrifugation at full speed will not affect yield or purity of the DNA. If, after centrifugation the sample has not completely passed through the membrane, centrifuge again until all the solution has passed through.

### 4. Protocol

1. Centrifuge the appropriate number of cells for 10 minutes at 1,000 x g. Avoid using more than  $5 \times 10^7$  mammalian cells per preparation.
2. Remove the supernatant and discard.
3. Resuspend the pellet in PBS to a final volume of 200 µL.
4. To each 200 µL sample of resuspended cells in PBS, add 20 µL of proteinase K.
5. *Optional RNase digestion steps: (If you do not wish to perform RNase digestion, move on to step 8)*
  - a. Add 4 µL RNase A (100 mg/mL) to the 220 µL sample of cells in PBS and Proteinase K. Pulse vortex to mix for 15 seconds.
  - b. Incubate for 2 minutes at room temperature.
  - c. Centrifuge for 5 seconds to remove drops from the inside of the lid.
6. Add 200 µL of Buffer AL and pulse vortex thoroughly for 15 seconds.
7. Incubate at 56 °C for 10 minutes to complete lysis.
8. Add 200 µL of 100% non-denatured ethanol to the sample and pulse vortex for 15 seconds to mix.
9. Centrifuge again for 5 seconds to remove drops from the lid.  
*(Thoroughly mix at each stage to obtain a homogenous solution).*
10. Apply the mixture, including any precipitate, to the NAB Nanosep device insert inside a receiver tube. If the sample volume is greater than 500 µL, apply half the sample to the NAB Nanosep device insert, perform steps 11 and 12 and then repeat with the second half of the sample.
11. Close the cap and centrifuge for 60 seconds at 10,000 – 14,000 x g.
12. Discard the flow-through and re-use the receiver tube for the next step.
13. Open the NAB Nanosep device and add 500 µL Buffer AW1.

14. Close the cap and centrifuge for 60 seconds at 10,000 – 14,000 x g.
15. Discard the collection tube and flow-through and transfer the NAB Nanosep device insert into a clean collection tube (provided).
16. Add 500 µL Buffer AW2 to the device insert.
17. Close the cap and centrifuge for 3 minutes at 10,000 – 14,000 x g.
18. Discard the flow-through and re-use the receiver tube for the next step.
19. Centrifuge for 60 seconds at 10,000 – 14,000 x g to remove possible Buffer AW2 carry-over.
20. Discard the receiver tube with the flow-through and transfer the device insert to the final, clean collection tube (provided).
21. To elute the DNA, add 100 µL of DNase-free distilled water (or Buffer AE) and incubate at room temperature for 5 minutes.
22. Centrifuge for 60 seconds at 10,000 – 14,000 x g.
23. *Optional: For maximum DNA yield, repeat the above elution steps with a further 100 µL of DNase-free distilled water.*

#### **Storage of DNA**

If you intend to store the purified DNA, we recommend elution with Buffer AE from Qiagen and storage at -30 °C to -15 °C. As high pH or EDTA can affect sensitive downstream applications, water can be used for elution. However, ensure that the pH of the water is a minimum of 7.0 as acid hydrolysis will occur to DNA stored in water which will result in degradation.

#### **Quantification of DNA**

DNA concentration can be determined by measuring the absorbance at 260 nm ( $A_{260}$ ) in a spectrophotometer (See details below). For small quantities of DNA however, it can be difficult to determine these amounts photometrically. Smaller quantities of DNA can be accurately quantified using fluorometric quantification.

#### **Spectrophotometric quantification of DNA**

$A_{260}$  readings should be greater than 0.10 and lower than 1.0 to ensure significance. An absorbance reading of 1.0 at 260 nm corresponds to 50 µg of DNA per mL. This is only valid for measurements at neutral pH however. As a result, if it is necessary to dilute the DNA sample, ensure that the dilution buffer is of neutral pH.




**Corporate Headquarters**  
25 Harbor Park Drive  
Port Washington, New York 11050

*Filtration. Separation. Solution.<sup>SM</sup>*

---

**Visit us on the Web at [www.pall.com/lab](http://www.pall.com/lab)**  
**E-mail us at [LabCustomerSupport@pall.com](mailto:LabCustomerSupport@pall.com)**

---

© 2020 Pall Corporation. Pall,  and Nanosep are trademarks of Pall Corporation. ® indicates a trademark registered in the USA. *Filtration. Separation. Solution.* is a service mark of Pall Corporation.