## Acrodisc® MS Syringe Filters

For LCMS Sample Preparation and Filtration of Organic or Aqueous Solvent-Based Samples

- Certified for Liquid Chromatography/Mass Spectrometry (LCMS)
- Extremely low levels of extractables
- Compatible with virtually all solvents
- Protects instruments against particulate build-up

### WWPTFE Membrane

The WWPTFE membrane is a water-wettable polytetrafluoroethylene membrane that is LCMS-certified. Using the Acrodisc MS syringe filter with WWPTFE membrane is ideal for LCMS sample prep. It offers extremely low extractable levels with good flow rates and retention efficiency. The Acrodisc MS syringe filter with WWPTFE membrane aids in obtaining accurate experimental results which saves time and money. Furthermore, the WWPTFE membrane provides excellent chemical compatibility with both organic and aqueous solvents.

### Ordering Information

#### Acrodisc MS Syringe Filters

Part Number	Description	Pkg
MS-3201	Acrodisc MS syringe filter, WWPTFE membrane, 0.2 µm, 25 mm	50/pkg

### Specifications

#### Materials of Construction

Filter Media: WWPTFE (water-wettable polytetrafluoroethylene) membrane  
Housing: HDPE (high density polyethylene)

**Pore Size**  
0.2 µm

**Effective Filtration Area**  
3.9 cm<sup>2</sup>

**Typical Sample Volume**  
< 150 mL

#### Inlet/Outlet Connections

Female threaded luer inlet, male slip luer outlet

#### Maximum Operating Pressure/Temperature

60 psi at 21-24 °C (70-75 °F)



30 psi at 55 °C (131 °F)

#### Typical Water Flow Rate

140 mL/min at 30 psi

### LCMS Certification

Pall Life Sciences certifies that all lots of the WWPTFE membrane used in the Acrodisc MS syringe filters have been tested according to established LCMS techniques to insure low levels of extractables.

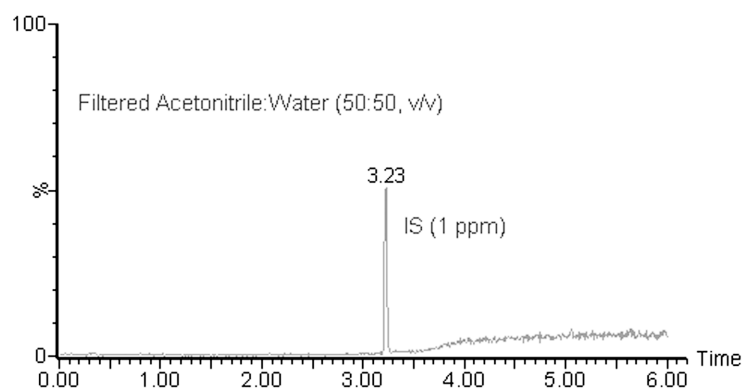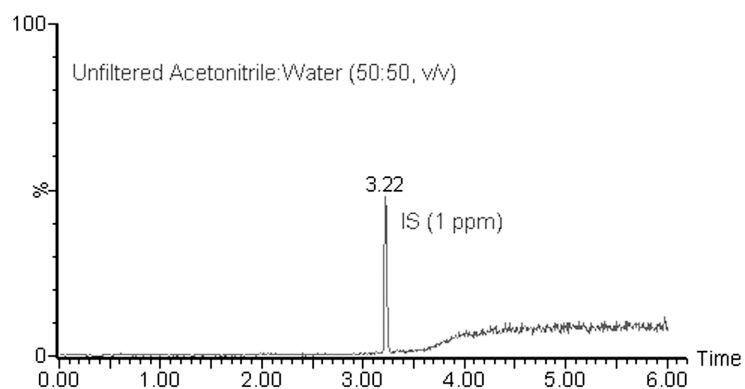### Test Parameters

Detection - Mass Spec.	ES+ Mode
	Mass Range: 100.00 - 1400.0
Column	Waters Acquity UPLC® BEH
	1.7 µm – 2.1 x 50 mm
Mobile Phase A	Methanol with 0.01% Formic Acid
Mobile Phase B	Water with 0.01% Formic Acid
Acquisition Program	Gradient - 6.0 minutes

#### IS: Internal Standard

The data provided above was obtained using the described conditions. Using other extraction methods may reduce or increase the extractable levels.

### Instructions for Use

Before filling the syringe with sample, draw approximately 1 mL of air into the syringe. This will allow the air to follow the sample out of the syringe. This "air purge" minimizes fluid retention within the filter device.

Fill a syringe with the solution to be filtered. To prevent accidental contamination, we recommend a glass syringe. CAUTION: Use of syringes smaller than 10 mL can generate excessive pressure on the filter, which may exceed the maximum operating pressure.

Holding the filter device in one hand and the filled syringe in the other, secure (without excessive force) the filled syringe to the filter device with a twisting motion.

Apply gentle pressure to begin filtration. (A gentle pressure helps ensure maximum throughput.) CAUTION: As the filter removes higher levels of particulates, filtration will become more difficult (the syringe plunger will be harder to move) and pressure will rapidly increase on the filter. Change filters when resistance becomes excessive.

**NOTE:** These filters are for **single use only**.

**Please avoid contact with potential sources of extractables during filtration and storage.**

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## WARNING

Employment of the products in applications not specified, or failure to follow all instructions contained in this product information insert, may result in improper functioning of the product, personal injury, or damage to property or the product. See Statement of Warranty in our most recent catalog.

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<b>Pall Life Sciences</b> 600 South Wagner Road Ann Arbor, MI 48103-9019 USA	<b>Australia</b> , Cheltenham, VIC, 03 9584 8100 <b>Austria</b> , Wien, 00 1 49 192 0 <b>Brazil</b> , São Paulo, SP, +55 11 5501 6000 <b>Canada</b> , Ontario, 905-542-0330 <b>Canada</b> , Québec, 514-332-7255 <b>China</b> , P.R., Beijing, 86-10-6780 2288 <b>France</b> , St. Germain-en-Laye, 01 30 61 32 32 <b>Germany</b> , Dreieich, 06103-307 333 <b>India</b> , Mumbai, 91 (0) 22 67995555 <b>Italy</b> , Buccinasco, +3902488870.1 <b>Japan</b> , Tokyo, 03-6901-5800 <b>Korea</b> , Seoul, 82-2-560-8711 <b>Malaysia</b> , Selangor, +60 3 5569 4892 <b>Poland</b> , Warszawa, 22 510 2100 <b>Russia</b> , Moscow, 5 01 787 76 14 <b>Singapore</b> , 65 6 389-6500 <b>South Africa</b> , Johannesburg, +27-11-2662300 <b>Spain</b> , Madrid, 91-657-9876 <b>Sweden</b> , Lund, (0)46 158400 <b>Switzerland</b> , Basel, 061-638 39 00 <b>Switzerland</b> , Fribourg, +41 (0)26 350 53 00 <b>Taiwan</b> , Taipei, 886 2 2545 5991 <b>Thailand</b> , Bangkok, 66 2937 1055 <b>United Kingdom</b> , Farlington, 02392 302600
1.800.521.1520 USA and Canada (+)800.PALL.LIFE Outside USA and Canada 1.734.665.0651 phone 1.734.913.6114 fax	
<b>Visit us on the Web at <a href="http://www.pall.com/lab">www.pall.com/lab</a></b>	
<b>E-mail us at <a href="mailto:LabSupport@pall.com">LabSupport@pall.com</a></b>	
© 2012, Pall Corporation. Pall,  , and Acrodisc are trademarks of Pall Corporation. ® indicates a trademark registered in the USA. ♦Waters Acquity UPLC is a trademark of Waters Corporation.	