MicroFunnel™ ST Sterility Testing Funnel

- Ideal for use in isolation chambers
- Disposable, ready-to-use filter funnel ideal for sterility testing applications
- Individually bagged funnels in overpack bag for easy transport and assured resistance to sterilizing agents used in isolation chambers
- Meets USP-24 requirements for Sterility Testing <71>

Ordering Information

Prod. No.	Description	Packaging
4851	MicroFunnel ST filter funnel with 0.2 µm Supor® membrane, gridded, 100 mL capacity	40/pkg*
4750	MicroFunnel ST filter funnel with 0.45 µm Supor membrane, plain, 100 mL capacity	40/pkg
4751	MicroFunnel 300ST filter funnel with 0.45 µm Supor membrane, plain, 300 mL capacity	20/pkg**
4811	MicroFunnel ST filter funnel with 0.45 µm GN-6 Metrical® membrane, gridded, 100 mL capacity	40/pkg
4812	MicroFunnel 300ST filter funnel with 0.45 µm GN-6 Metrical membrane, gridded, 300 mL capacity	20/pkg

*100 mL capacity funnels packaged 10 individually bagged funnels per overpack bag, 4 overpack bags per box

**300 mL capacity funnels packaged 5 individually bagged funnels per overpack bag, 4 overpack bags per box

Membranes

GN-6 Metrical membrane: A 0.45 µm mixed cellulose ester membrane ideal for use with aqueous solutions

Supor membrane: 0.2 or 0.45 µm polyethersulfone membrane ideal for low binding of antibiotics or other inhibitory substances

Specifications

Materials of Construction

Filter Media:

PN 4750, 4751, 4851: Supor (polyethersulfone) membrane

PN 4811, 4812: GN-6 Metrical (mixed cellulose ester) membrane

Support Pad: Cellulose

Funnel and Base: Polypropylene

Funnel Cover: Polystyrene

Bag and Overpack Bag: Polyethylene

Adapter: Polyethylene


Effective Filtration Area

13.46 cm²

Maximum Vacuum

635 mm Hg (25 in. Hg) (Vacuum use only)

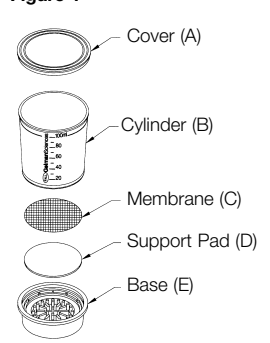Figure 1

Specifications (cont.)

Sterilization

Sterilized by gamma irradiation

Dimensions

100 mL Funnels

Height: 7.62 cm (3.00 in.) with cover
7.32 cm (2.88 in.) without cover

Diameter: 6.05 cm (2.38 in.) with cover
5.74 cm (2.26 in.) without cover

300 mL Funnels

Height: 9.09 cm (3.58 in.) with cover
8.94 cm (3.52 in.) without cover

Diameter: 8.84 cm (3.48 in.) with cover
8.53 cm (3.36 in.) without cover

Instructions for Use

Please consult the appropriate compendial methods for information and guidelines on selection and validation of rinse fluid.

Transfer of test materials into isolator chamber:

1. Place overpack bag containing individually bagged funnels into transfer chamber.
2. Run sterilization cycle.
3. Using sterilized scissors, cut open the overpack bag being careful to avoid damage to the individual bags contained within.
4. Transfer individually bagged product into testing chamber.

Process the sample:

5. Remove the sterile, disposable funnel by tearing bag open at the notch and place funnel onto vacuum manifold set-up as illustrated in Figure 2.
6. Remove cover without touching the inner surface and pour sample into the funnel. Replace cover.
7. Turn on the vacuum and draw the sample through the filter.
8. Remove the cover and rinse funnel thoroughly with a volume of sterile, buffered solution appropriate to your sample. Replace cover until rinse solution is drawn completely through the filter.
9. Separate the cylinder from the base by removing cover and gently squeezing near the top of the cylinder.
10. Carefully remove the membrane filter from the base with smooth-tipped forceps.
11. Either submerge the entire membrane in nutrient broth medium (Fluid Thioglycollate or Fluid Soybean Casein Digest medium) or using sterile scissors, cut the membrane into two equal parts and submerge each half in separate nutrient broth media.
12. Incubate nutrient broth media as per pharmacopeial specifications.

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
Figure 2

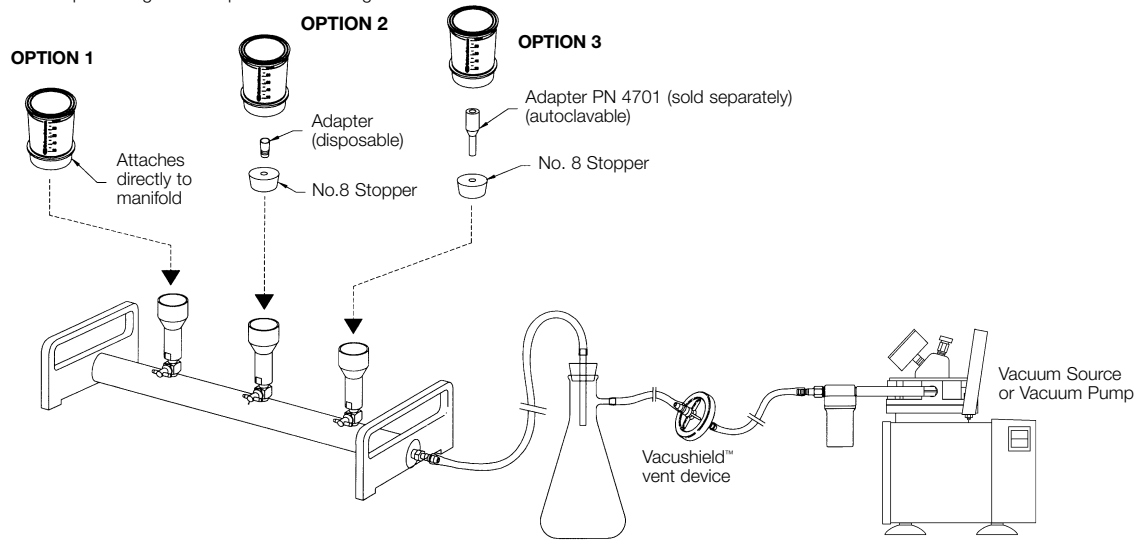Filtration manifold setup detailing various options of attaching filter units to manifold.

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Complementary Products		
Prod No.	Description	Packaging
Hardware and Accessories		
15402	3-Port Aluminum Manifold	1/pkg
15403	6-Port Aluminum Manifold	1/pkg
13157	Vacuum/Pressure Pump, 115 V	1/pkg
13158	Vacuum/Pressure Pump, 230 V CE	1/pkg
13160	Vacuum Pump, 115 V	1/pkg
51147	Stainless Steel Forceps	1/pkg
4701	Autoclavable Funnel Adapter	3/pkg
82728	No. 8 Rubber Stopper	1/pkg
4402	Vacushield™ Vent Device	3/pkg
4713	Vacuum Pump Funnel Adapter	1/pkg
Additional MicroFunnel™ Filter Funnels		
4800	MicroFunnel filter unit with 0.45 µm GN-6 Metricel®, gridded, 100 mL capacity, individually bagged	50/pkg
4801	MicroFunnel NB filter unit with 0.45 µm GN-6 Metricel, gridded, 100 mL capacity	50/pkg
4803	MicroFunnel SP filter unit with 0.2 µm Supor 200®, gridded, 100 mL capacity, individually bagged	50/pkg
4804	MicroFunnel filter unit with 0.45 µm GN-6 Metricel, gridded, 100 mL capacity, individually bagged	200/pkg
4805	MicroFunnel YMB filter unit with 0.45 µm Metricel Black, gridded, 100 mL capacity, individually bagged	50/pkg
4806	MicroFunnel SNG filter unit with 0.2 µm Supor 200, plain, 100 mL capacity, individually bagged	50/pkg
4810	MicroFunnel LP filter unit with 0.45 µm GN-6 Metricel, gridded, 100 mL capacity, individually bagged	50/pkg
4815	MicroFunnel 300 filter unit with 0.45 µm GN-6 Metricel, gridded, 300 mL capacity, individually bagged	20/pkg
4817	MicroFunnel 300 YMB filter unit with 0.45 µm Metricel Black, gridded, 300 mL capacity, individually bagged	20/pkg
4818	MicroFunnel 300 SP filter unit with 0.2 µm Supor 200, gridded, 300 mL capacity, individually bagged	20/pkg
Additional MicroFunnel Plus Filter Funnels		
4807	MicroFunnel Plus filter funnel with 0.45 µm GN-6 Metricel membrane, 100 mL capacity, individually bagged	50/pkg
4808	MicroFunnel Plus filter funnel with 0.45 µm Metricel Black membrane, 100 mL capacity, individually bagged	50/pkg
4809	MicroFunnel Plus filter funnel with 0.2 µm Supor membrane, 100 mL capacity, individually bagged	50/pkg

WARNING

Employment of the products in applications not specified, or failure to follow all instructions contained in this product information insert, may result in improper functioning of the product, personal injury, or damage to property or the product. See Statement of Warranty in our most recent catalog.

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Pall Life Sciences

600 South Wagner Road
Ann Arbor, MI 48103-9019 USA

For ordering or technical information:

In USA and Canada
Tel: 734-665-0651
800-521-1520
Fax: 734-913-6114
Outside USA and Canada
+800 PALL LIFE
+800 7255 5433

Visit Pall Life Sciences on the Web at www.pall.com/Lab or e-mail us at Lab@pall.com

Offices:
Australia, Cheltenham, VIC, 03 8586 8150
Austria, Wien, 00 1 49 192 0
Canada, Ontario, 905-542-0330
Canada, Québec, 514-332-7255
China, P.R., Beijing, 86-10-8458 4010
France, St. Germain-en-Laye, 01 30 61 39 92
Germany, Dreieich, 06103-307 333
India, Mumbai, 91 (0) 22 55995555
Italy, Milano, 02 47 79 61
Japan, Tokyo, 03-6901-5800
Korea, Seoul, 82-2-560-7834
Malaysia, Selangor, +60 3 5569 4892
New Zealand, Hamilton, +64 7 957 9510
Poland, Warszawa, 22 510 2100
Russia, Moscow, 5 01 787 76 14
Singapore, (65) 389-6500
South Africa, Johannesburg, +27-11-2662300
Spain, Madrid, 91-657-9876
Sweden, Lund, (0)46 158400
Switzerland, Basel, 061-638 39 00
Taiwan, Taipei, 886 2 2545 5991
Thailand, Bangkok, 66 2937 1055
United Kingdom, Farington, 02392 302600