



O-Ring Replacements for Aluminum Manifold Valves

- Kit contains enough o-rings to service a 3-place manifold

Ordering Information

Prod. No.	Description	Packaging
88160	O-ring replacement kit (3 size -060, 6 size -010)	1 kit

Replacement Instructions

1. Carefully remove the c-clip by evenly pushing the two ends of the c-clip until it slides off perpendicular to the end of the valve stem.
2. Pull the stem out of the valve case. Note the orientation of the flat side on the valve stem in relationship to the stop on the valve case.
3. Remove old o-rings and clean valve of any residue if necessary.
4. Replace with new o-rings: one (1) size -060 and two (2) size -010 per valve. Silicone lubricant will aid in getting the new o-rings onto the valve stem and enable the assembled valve to move smoothly.
5. Carefully press the valve stem into the valve case with the flat side on the valve stem facing the stop on the valve case. Work the stem gently back and forth as you press it into place so the o-rings don't get pinched and damaged.
6. Slide c-clip back into place on the end of the valve stem.

Complementary Products

Prod. No.	Description	Packaging
15402	3-place Aluminum Manifold	1/pkg
15403	6-place Aluminum Manifold	1/pkg

WARNING

Employment of the products in applications not specified, or failure to follow all instructions contained in this product information insert, may result in improper functioning of the product, personal injury, or damage to property or the product. See Statement of Warranty in our most recent catalog.

**PALL**[®]

Life Sciences

Pall Life Sciences

600 South Wagner Road
Ann Arbor, MI 48103-9019 USA

For ordering or technical information:

Tel: 734-665-0651

800-521-1520 (in USA)

Fax: 734-913-6114

**Visit Pall Life Sciences on the
Internet: www.pall.com/Lab or
e-mail us at: Lab@pall.com**

Offices:

Australia, Lane Cove, NSW, 02-9428 2333

Austria, Wien, 043-1-49 192 0

Canada, Ontario, 905-542-0330

Canada, Québec, 514-332-7255

China, P.R., Beijing, 86-10-8458 4010

France, St. Germain-en-Laye, 01 30 61 39 92

Germany, Dreieich, 06103-307 333

India, Mumbai, 91-22-5956050

Italy, Milano, 02-47-79-61

Japan, Tokyo, 3-3495-8319

Korea, Seoul, 2-569-9161

Poland, Warszawa, 22-835 83 83

Russia, Moscow, 095 787-76-14

Singapore, 65-389-6500

Spain, Madrid, 91-657-9876

Sweden, Lund, 46-0-46 15 84 00

Switzerland, Basel, 061-638 39 00

Taiwan, Taipei, 2-2545-5991

United Kingdom, Portsmouth, 023 92 302600

Pall, and , are trademarks of Pall Corporation. © indicates a registered trademark in the USA.