**TCLP Glass Fiber Acrodisc® Syringe Filter**

- Fast flow rates and high particulate holding capacities for prefiltration applications
- 0.7 µm, nominal rating is ideal for water and wastewater analysis

Ordering Information

Prod. No.	Description	Packaging
4495	TCLP Glass Fiber Acrodisc Syringe Filter, 0.7 µm, 25 mm	1000/pkg

Specifications**Materials of Construction**

Filter Media: TCLP Glass Fiber; binder free borosilicate glass fiber
Housing: Polypropylene

Pore Size

0.7 µm, nominal

Effective Filtration Area

2.8 cm²

Sample Volume

< 100 mL

Inlet/Outlet Connections

Ferrule Luer-Lok® inlet, standard male slip luer outlet

Typical Hold-up Volume (with Air Purge)

< 125 µL

Maximum Operating Temperature

55 °C (131 °F)

Maximum Operating Pressure

5.2 bar (75 psf)

Instructions for Use**English**

Before filling the syringe with sample, draw approximately 1 mL of air into the syringe. This will allow the air to follow the sample out of the syringe. This "air purge" minimizes fluid retention within the filter device.

Fill the syringe with the solution to be filtered. CAUTION: Use of syringes smaller than 10 mL can generate excessive pressure on the filter, which may exceed maximum operating pressure.

Hold the filter device in one hand and the filled syringe in the other, secure (without excessive force) the filled syringe to the filter device with a twisting motion. Apply gentle pressure to begin filtration. (A gentle pressure helps assure maximum throughput.) CAUTION: As the filter removes particulates, filtration will become more difficult (the syringe plunger will be harder to use) and pressure will rapidly increase on the filter. Change filters when resistance becomes excessive. Failure to change filter may result in housing rupture, which results in particulate contaminating the filtrate.

These filters are for **SINGLE USE ONLY**.

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